



Clean Water. Fertile Soil. Serious Science.

334 Hecla Street, Lake Linden, MI 49945 USA

Toll-free: 888/NITRATE

Int'l phone: 1.906.296.1115

Fax: 1.906.296.8003

Website: www.nitrate.com

Tech Line: 1.906.296.1130

Support: tech@nitrate.com

Data Sheet for Phosphate Test Kits

Phosphate reaction cuvettes do not produce a visible color change and must be read photometrically. Refer to the photometer manual if using the NECi Handheld photometer.

For users utilizing their own photometer, read the reactions at an absorbance of 360 nm. Below are typical curves for the phosphate test kits. Create a standard curve and plot samples against the curve to get phosphate results.

<i>Sample Number</i>	<i>Sample Description</i>	<i>ppm PO4/ Abs 360</i>
Example	Northwest corner of pond, collected 8/25/15	Close to, but less than 10ppm
1		
2		
3		
4		
5		

